MCE USA

Tel: 609-228-6898

Email: tech@MedChemExpress.com



Phalloidin dye conjugates

1 Contents

Cat. No.	Product Name	Ex (nm)	Em (nm)	Size	Package
HY-K0901	AMCA Phalloidin	344	432	300 T	30 μL
HY-K0902	Fluorescein Phalloidin	497	516	300 T	30 μL
HY-K0903	Rhodamine Phalloidin	552	578	300 T	30 μL

2 General Information

Phalloidin is a mushroom-derived toxin which can be used to specifically label F-actins (actin filaments) of the cytoskeleton. It binds to all variants of actin filaments in many different species of animals and plants. Typically, Phalloidin is used conjugated to a fluorescent dye. Phalloidin binds F-actins with high selectivity while fluorescent dye provides stable and bright fluorescence.

MCE offers three Phalloidin dye conjugates, AMCA Phalloidin (HY-K0901, blue fluorescence), Fluorescein Phalloidin (HY-K0902, green fluorescence) and Rhodamine Phalloidin (HY-K0903, orange fluorescence). These Phalloidin dye conjugates are 1000× stock solution in DMSO which is convenient for labeling, identifying and quantitating F-actins in in formaldehyde-fixed and permeabilized tissue sections, cell cultures or cell-free experiments.

3 Protocol

1. Prepare 1× Phalloidin dye conjugate working solution:

Add 1 µL of 1000× Phalloidin dye conjugate DMSO solution to 1 mL of PBS containing with 1% BSA.

Note 1: Different cell types might be stained differently. The concentration of Phalloidin dye conjugate working solution should be prepared accordingly.

Note 2: 1% bovine serum albumin (BSA) is used to reduce nonspecific background staining.

- 2. Stain the cells:
- 2.1 Wash cells 2-3 times with PBS. Fix the cells in 3.7% methanol-free formaldehyde solution in PBS for 10-30 minutes at room temperature.

Note: Avoid any methanol containing fixatives since methanol can disrupt actin during the fixation process. The preferred fixative is methanol-free formaldehyde.

- 2.2 Wash the fixed cells 2-3 times in PBS.
- 2.3 Optional: Add 0.1% Triton X-100 in PBS into fixed cells for 3–5 minutes to increase permeability. Wash the cells 2–3 times in PBS.
- 2.4 Add 100 µL/well (96-well plate) of Phalloidin dye conjugate working solution into the fixed cells, and stain the cells at room temperature for 20 to 90 minutes.
- 2.5 Rinse cells gently with PBS 2-3 times to remove excess Phalloidin dye conjugate.
- 2.6 Run fluorescence microscope or other equipments at desired Ex/Em wavelengths (AMCA Phalloidin, Ex/Em=344/432 nm; Fluorescein Phalloidin, Ex/Em=497/516nm; Rhodamine Phalloidin, Ex/Em=552/578 nm).

4 Storage

The solutions should be stable for at least 6 months if store at -20°C.

Protect the Phalloidin dye conjugates from light, and avoid freeze/thaw cycles.

5 Precautions

- 1. One Test(T) of Phalloidin dye conjugates is equivalent to 0.1 μ L of the DMSO stock solution (100 μ L working solution).
- 2. Phalloidin is toxic although the amount of toxin present in a vial could be lethal only to a mosquito (LD_{50} of phalloidin = 2 mg/kg). Please handle with care.
- 3. This product is for R&D use only, not for drug, household, or other uses.